

Blades Biological Ltd

COWDEN . EDENBRIDGE . KENT . UK . TN8 7DX

Telephone: (01342) 850242 . Fax: (01342) 850924

Email: info@blades-bio.co.uk Online: <http://www.blades-bio.co.uk/>

Brine Shrimp Eggs - *Artemia*

Brine shrimp eggs should be hatched into a container with a large surface area.

Fill the container to within 3cm of the top with **Marine Salt Solution** (see our current catalogue).

Sprinkle shrimp eggs evenly onto the water surface. For best results, use 1g of eggs to 1 litre of solution.

At room temperature, (21°C) the eggs should hatch in 24 to 36 hours.

To remove the shrimp from the solution, either strain them through a fine mesh or, using a source of light, attract them towards one spot and remove them using a pipette.

For feeding purposes, if necessary, rinse shrimp several times with fresh bottled water to remove all traces of salt.