

Blades Biological Ltd

COWDEN . EDENBRIDGE . KENT . UK . TN8 7DX

Telephone: (01342) 850242 . Fax: (01342) 850924

Email: info@blades-bio.co.uk Online: <http://www.blades-bio.co.uk/>

VINEGAR EEL CARE NOTES. *Anguilla Aceti*.

On receipt of your vinegar eel culture loosen the jar lid and place in an area out of direct sunlight at room temperature.

To subculture, pour half of the culture into two clean jars and fill with cider vinegar.

The cultures should be kept at 20'C to 30'C and fresh vinegar should be added to balance evaporation. The population of worms should become dense within a few weeks and the culture should then be sub cultured again.